# CATHEPSIN K

ENZYME IMMUNOASSAY FOR THE QUANTITATIVE DETERMINATION OF CATHEPSIN K  
IN SERUM OR CELL CULTURE SUPERNATANTS  
CAT. NO. BI-20432 12 X 8 TESTS

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For Research Purposes Only

FOR RESEARCH USE ONLY  
NOT FOR USE IN DIAGNOSTIC PROCEDURES

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**[www.bmgrp.com](http://www.bmgrp.com)**

## 1) INTRODUCTION

Cathepsin K is a highly expressed cysteine protease, secreted by osteoclasts. It is produced as a 329 AA pre-cursor, which is cleaved to its active form with a length of 215 AA (MW 23,495) by osteoclasts, when they are resorbing bone. Cathepsin K degrades the collagen matrix of the bone after removal of its mineral components. It is actually the most abundantly synthesised protein of the active, resorbing osteoclast and therefore the most specific marker of resorption activity. Cathepsin K plays a key role in tissue destruction, remodelling and cartilage breakdown of the bone. The high homology of Cathepsin K across various species (mouse 86%, rat 88%, pig 97%, rabbit 96%) should allow the use of this assay for studies in animal models for following diseases

### RESEARCH AREAS

- Primary and secondary osteoporoses
- Inflammatory processes which affect multiple regions of the skeleton
- Antiresorptive drug monitoring
- Bone turnover research

## 2) CONTENTS OF THE KIT

CONT	KIT COMPONENTS	QUANTITY
PLATE	Polyclonal sheep anti-Cathepsin K antibody coated microtiterstrips in strip holder, packed in alu bag with desiccant	12 x 8 tests
WASHBUF	Washing buffer 20x concentrate, natural cap	1 x 50 ml
ASYBUF	Assay Buffer, red cap, ready to use	1 x 25 ml
STD	Standards, (0; 11; 33; 100; 300 pmol/l), white caps, lyophilised with blue dye	5 x lyophilised
CTRL	Control, yellow cap, lyophilised, concentration after reconstitution see label	1 x lyophilised
CONJ	Conjugate, (polyclonal anti-Cathepsin K-HRP), amber cap, ready to use	1 x 22 ml
SUB	Substrate (TMB solution), blue cap, ready to use	1 x 22 ml
STOP	Stop solution, white cap, ready to use	1 x 7 ml

## 3) ADDITIONAL MATERIAL ADDED TO THE KIT

- 1 self-adhesive aluminium plastic film
- QC protocol
- Protocol sheet
- Instruction manual for use

## 4) EQUIPMENT REQUIRED BUT NOT SUPPLIED

- Precision pipettes calibrated to deliver 50, 200 and 1000 µl, disposable tips
- ELISA reader for absorbance at 450 nm (Reference 620 nm)
- Graph paper or software for calculation of results
- Plate washer is recommended for washing
- Distilled or deionised water

## 5) REAGENTS AND SAMPLE PREPARATION

Sample collection:

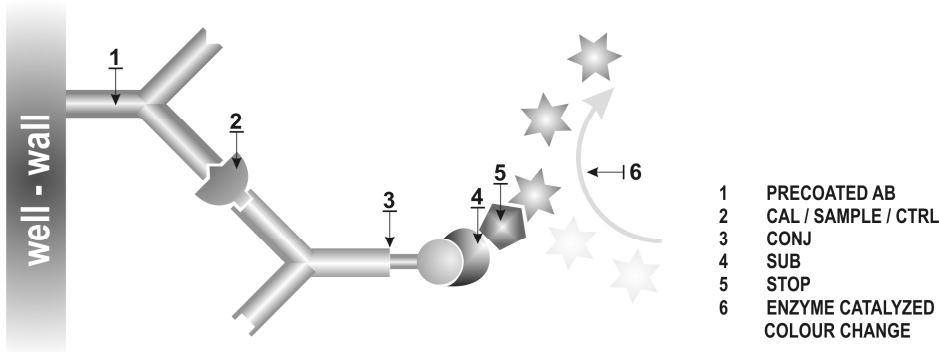
Freshly collected blood samples must be centrifuged and tested within 90 minutes if handled at room temperature. If handled at +4°C Cathepsin K is stable up to 2 days. For long term storage keep the samples at least at -20°C or better at -70°C. Up to 4 freeze/thaw cycles do not change the results. Lipemic or hemolyzed samples may give erroneous results. Samples should be mixed well before assaying.

Cell culture supernatants do not need any special pre-treatment. They can also be read against the supplied STD (standards).

#### Reconstitution / Handling:

- WASHBUF (Wash buffer): Dilute the concentrate 1:20 (1+19) e.g. 50 ml WASHBUF + 950 ml distilled water. Crystals in the buffer concentrate will dissolve at room temperature. The diluted WASHBUF is stable at 2-8°C until expiry date stated on label. Use only diluted WASHBUF (Wash buffer) for the assay performance.
- STD (Standard) and CTRL (Control): Pipette 1ml of ASYBUF (Assay buffer) to the vial. Leave at room temperature (18-26°C) for 20 min. **MIX WELL** on a Vortex Mixer. Aliquot the standards into Eppendorf tubes and freeze the standards as soon as possible after usage. Do not let them stand at room temperature. Reconstituted standard and control are stable at -20°C or -70°C until expiry date stated on the label.  
Avoid freeze – thaw cycles.

### 6) PRINCIPLE OF THE ASSAY

### 7) ASSAY PROTOCOL

All reagents and samples must be at room temperature (18-26°C) before use in the assay.
Mark position for BLANK/STD/SAMPLE/CTRL (Blank/Standard/Sample/Control) on the protocol sheet.
Take microtiter strips out of the alu bag, take a minimum of one well as Blank. Store unused strips with desiccant at 2-8°C in the alu bag. Strips are stable until expiry date stated on the label.
Add 50 µl STD/SAMPLE/CTRL (Standard/Sample/Control) in duplicate into respective wells, except blank.
Add 200 µl CONJ (anti Cathepsin K-HRPO) into each well, except blank, swirl gently.
<b>Cover tightly and incubate over night (20-24h) at room temperature (18-26°C) in the dark.</b>
Aspirate and wash wells 5x with 300 µl diluted WASHBUF (Wash buffer), remove remaining WASHBUF by hitting plate against paper towel after the latest wash.
Add 200 µl SUB (Substrate) into each well.
<b>Incubate for 30 min at room temperature (18-26°C) in the dark.</b>
Add 50 µl STOP (Stop solution) into each well.
Measure absorbance immediately at 450 nm with reference 620 nm, if available.

### 8) CALCULATION OF RESULTS

Use commercially available software or graph paper for calculation of the results. Subtract the blank extinction from all values and construct the standard curve from the standard values. Obtain the concentration of the samples from this standard curve. Respective dilution factors have to be considered.

The quality control protocol supplied with the kit shows the results of the final release QC for each kit. Data for optical density obtained by customers may differ due to various influences and/or due to the normal decrease of signal intensity during shelf life. However, this does not affect validity of results as long as an optical density of 1.50 is obtained for the standard with the highest concentration.

### 9) ASSAY CHARACTERISTICS

Normal range:	Median: 8.7 pmol/l (n = 30) It is recommended to establish the normal range for each laboratory.
Standard range:	0 to 300 pmol/l
Sample volume:	50 µl human serum or cell culture supernatant
Detection Limit: (0 pmol/l + 3 SD):	1.1 pmol/l
Incubation time:	Over night (20-24h) / 30 min

### 10) PRECISION

Intra-Assay (n=16)			Inter-Assay (n=10)		
Mean (pmol/l)	72	298	Mean (pmol/l)	62	276
SD	3	18	SD	4	22
CV%	4%	6%	CV%	6%	8%

### 11) TECHNICAL HINTS

- Do not mix or substitute reagents with those from other lots or sources.
- Do not mix stoppers and caps from different reagents or use reagents between lots.
- Do not use reagents beyond expiration date. Protect reagents from direct sunlight.
- Substrate solution should remain colourless until added to the plate.
- To ensure accurate results, proper adhesion of plate sealers during incubation steps is necessary.
- Avoid foaming when mixing reagents.

### 12) PRECAUTIONS

All test components of human source were tested with 3<sup>rd</sup> generation tests against HIV Ab and HBsAg; and were found negative. Nevertheless, they should be handled and disposed as if they were infectious.

All liquid reagents contain 0.01% Proclin 300 as preservative.

Proclin 300 is not toxic in concentrations used in this kit. It may cause allergic skin reactions – avoid contact with skin or eyes.

- Do not pipette by mouth.
- Do not eat, drink, smoke or apply cosmetics where reagents are used.
- Avoid all contact with the reagents by using gloves. Sulfuric acid is irritating to eyes and skin. Flush with water if contact occurs. Avoid contact with skin and mucous. Irritations are possible – Flush with water after contact!!

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# **BI-20432 CATHEPSIN K**

## **ASSAY PROTOCOL AND CHECKLIST**

- Bring all reagents to room temperature (18-26°C).
- Prepare reagents and samples as instructed.
- Bring unused and prepared components to the storage temperature mentioned in the package insert.
- Take microtiter strips out of the foil bag and mark positions on the protocol sheet.
- Add 50 µl STD/SAMPLE/CTRL (standard/ sample/control) into all wells, except blank.
- Add 200 µl CONJ (Conjugate) into each well, except blank, swirl gently.
- Cover tightly and incubate over night (20-24h) at room temperature (18-26°C) in the dark.**
- Aspirate and wash wells with 300 µl diluted WASHBUF (Wash buffer) five times. Remove remaining buffer by hitting plate against paper towel.
- Add 200 µl SUB (Substrate) into each well.
- Incubate for 30 minutes at room temperature (18-26°C), in the dark.**
- Add 50 µl STOP (Stop solution) into each well.
- Read Optical Density at 450 nm with reference 620 nm, if available.